
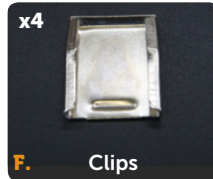


TrippNT Core CL™ Cart

Assembly Guide (v.3.1)

Thank you for choosing TrippNT for your storage and organization needs. Please check to make sure you received your cart in good shape and all the necessary components are included. This guide will show you step by step how to assemble your Core CL™ Cart.


Included Components

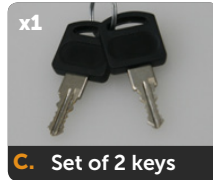


If you need any assistance please call TrippNT at 1.800.874.7768 or email us at Customerservice@TrippNT.com.

TrippNT Core CL™ Cart

Unpacking the Cart

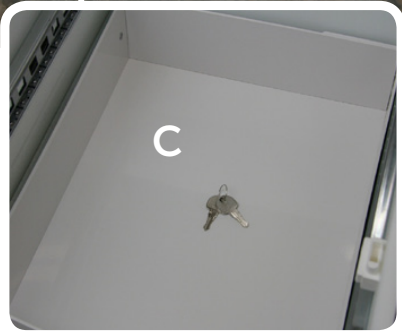


Locate the keys (C) taped to the second drawer.

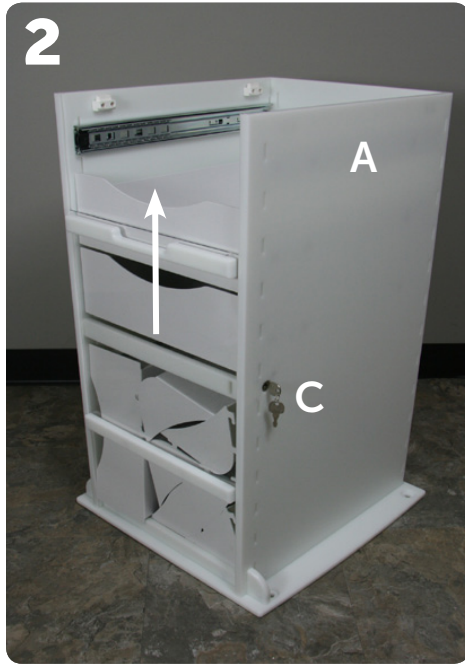

Unlock sliding door with keys (C).
Slide door up.



Unpack components and inspect with the included component list to ensure all items have been included.

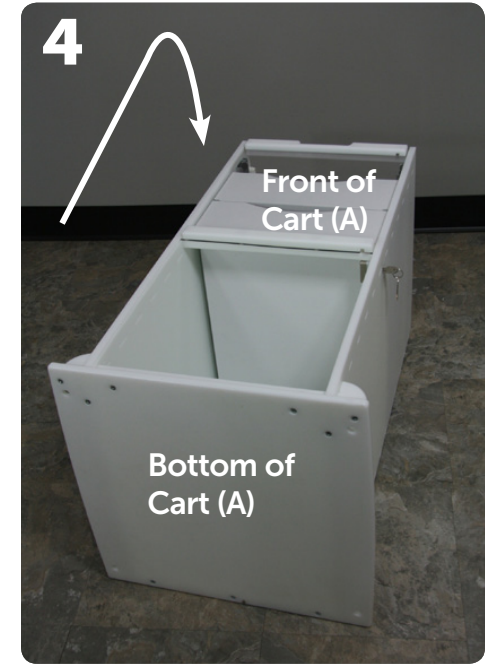

Gently lay cart body (A) on its back with the drawers and door facing up towards the ceiling.

Care Instructions for Your Cart

This product can be cleaned with a solution of water or with a mild organic solvent. Use of strong solvents is not recommended.

Accessories and Replacement Parts


A variety of accessories and replacement parts are available at:

[TrippNT.com/CartAccessories](https://www.trippnt.com/CartAccessories) or from your normal TrippNT Distributor

TrippNT Core CL™ Cart


Installing the Casters

5

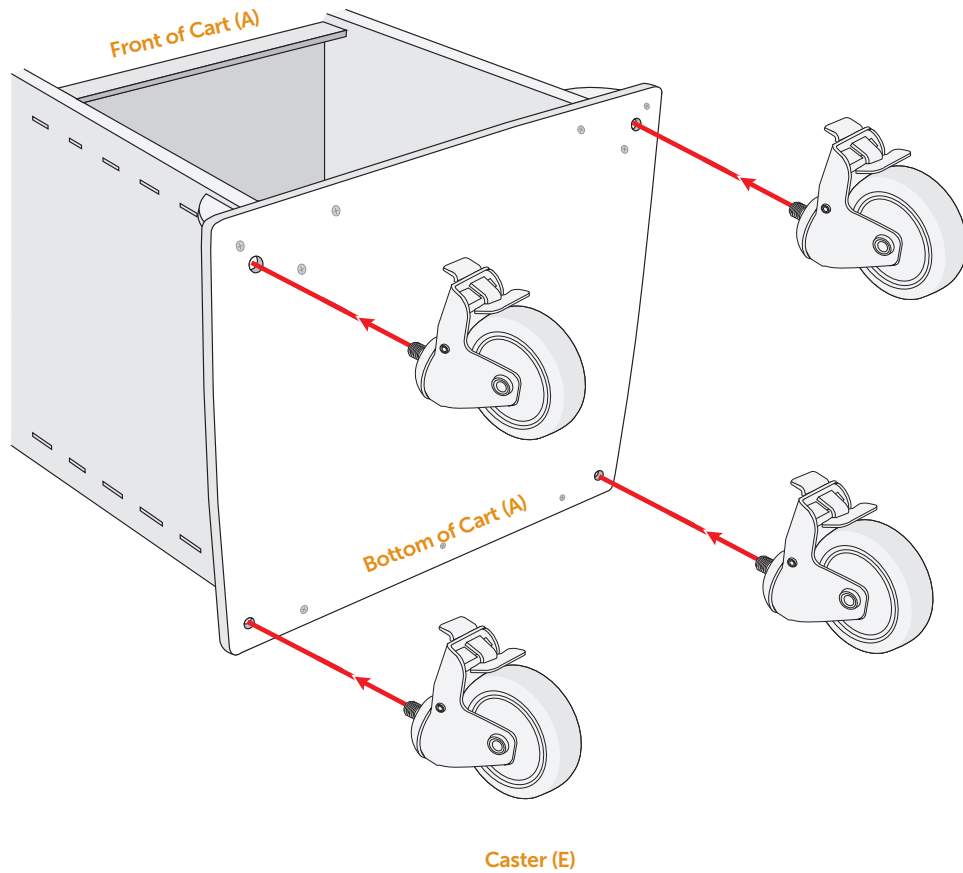

IMPORTANT CASTER INSTALLATION INSTRUCTIONS

Castors must be assembled correctly. The caster will stay on tight if installed correctly.

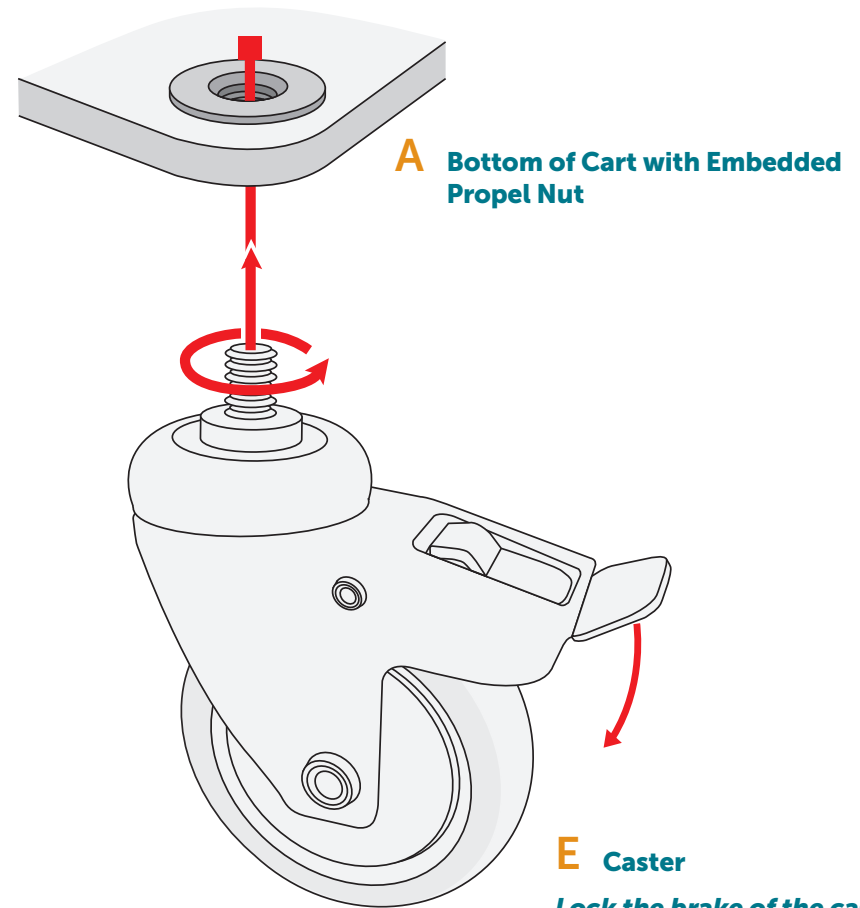

Lock the brake of the caster. Turn the caster in a clockwise direction until it is snug.



Keep the brake on all of the casters locked until assembly is complete.

TrippNT Core CL™ Cart


Assembling the Cart




Return cart to upright position. Ensure the lock brake on the front casters is engaged to prevent movement while assembling.

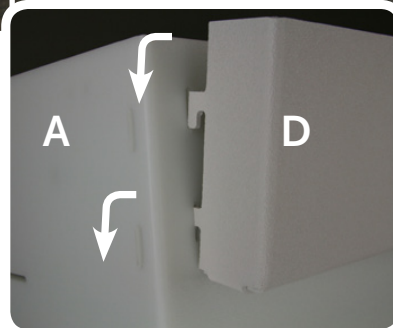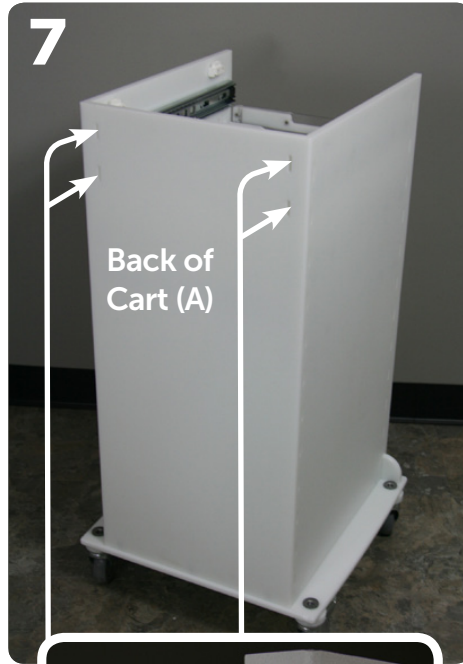

On the back of the cart body (A) Insert the tabs of tool holder (D) into the four slots and press down to secure.

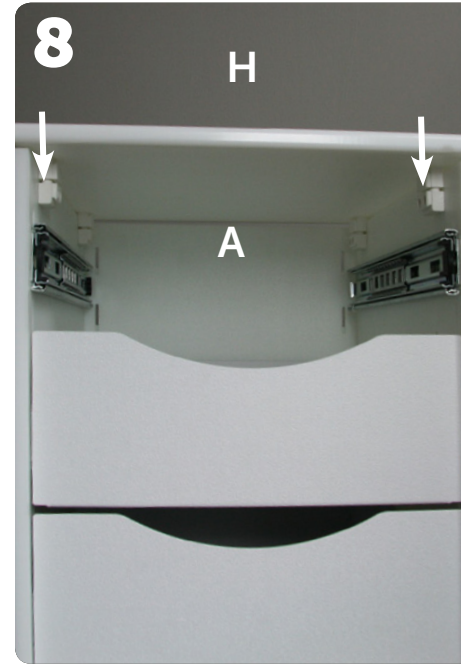


Install the cart top (H) using the plastic quick fittings. The fittings will snap when the cart top is in place. **Please note the side of the cart top with the groove for the door should face towards the front of the cart.**

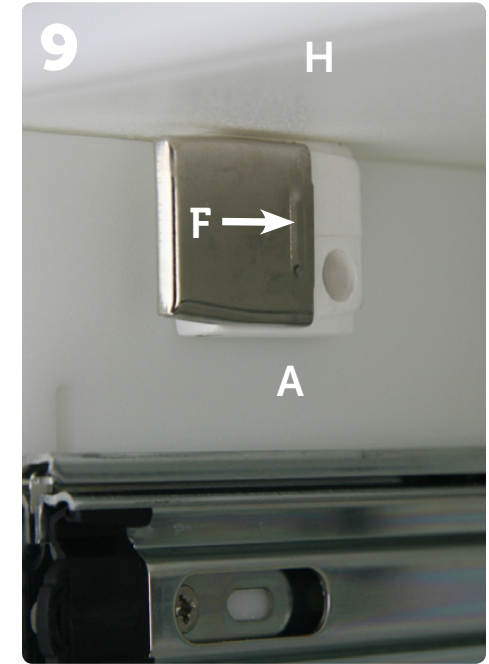



Slide metal clips (F) into position to secure the cart top (H) to the cart body (A).

Note:

Install the tool holder before installing the top of the cart.

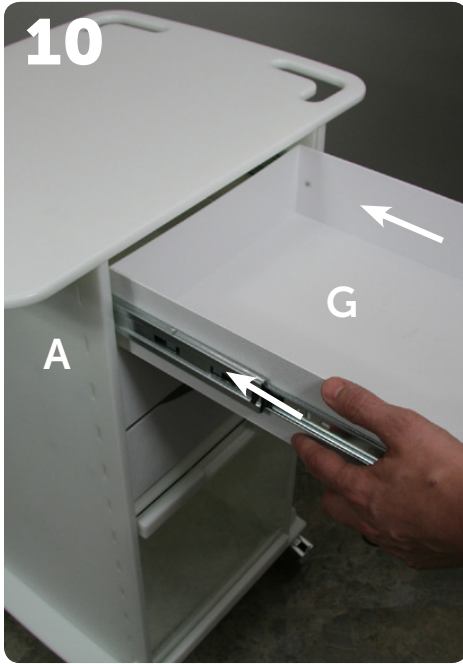

Align the tracks of the top drawer (G) with the tracks on the inside of the cart body. Press the top drawer completely in to the cart to snap it into place.

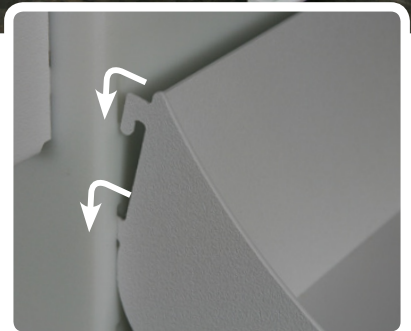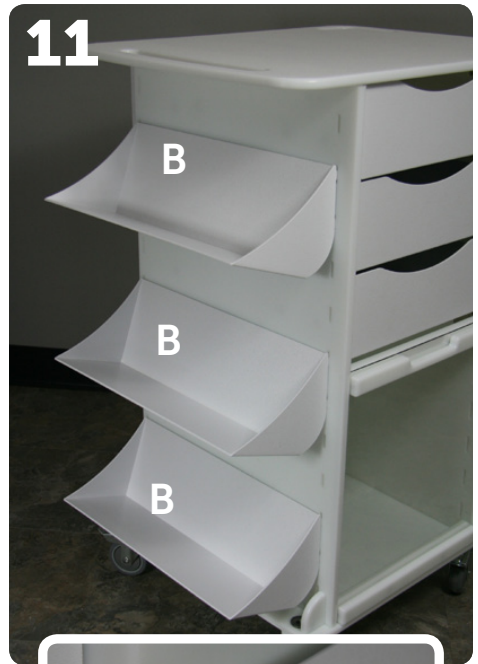

Insert the tabs from the side shelves (B) into the slots as shown at your desired height. Repeat on the other side and your cart is complete.